

Dec 30, 2017

DNA Analysis by Native Polyacrylamide Gel Electrophoresis and Infrared Fluorescence Imaging V.2

Michael Van Dyke¹, James Cox¹

¹Department of Chemistry and Biochemistry, Kennesaw State University

Working dx.doi.org/10.17504/protocols.io.mcy2xw



Michael Van Dyke
Department of Chemistry and Biochemistry, Kennesaw State Uni...



ABSTRACT

Short DNA probes, for example, those used in characterizing protein-DNA complexes, have historically been radiolabeled with ³²P to allow their detection and quantitation following native polyacrylamide gel electrophoresis (PAGE). For reasons of economy or safety, alternative means of DNA detection need to be used. Described here is the resolution of 5' IRD700 and IRD800 end-labeled DNAs by native PAGE and their visualization and quantitation by IR fluorescence imaging.

GUIDELINES


Protocol development supported by grants from the NIH ([1R15GM104833-01](#)) and NSF ([MCB-1714778](#)).

SAFETY WARNINGS


Acrylamide dust is an inhalation hazard and cumulative neurotoxin. Best to wear a protective face mask when weighing.

Handcast native polyacrylamide gels

- 1 We routinely use the Bio-Rad Mini-PROTEAN® vertical electrophoresis system to cast and run 10-well 1.0 mm-thick 9.5%:0.5% polyacrylamide:bisacrylamide minigels in ½X TBE buffer [44.5 mM Tris base, 44.5 mM boric acid, 1 mM EDTA]. Needed materials: Bio-Rad casting stand, two casting frames, two short plates, two 1.0 mm spacer plates, two 1.0 mm 10-well combs, and one Mini-PROTEAN® vertical electrophoresis cell with buffer dam. Needed reagents: 5x TBE buffer, 38:2 acrylamide:bisacrylamide, 10% ammonium persulfate (APS), and tetramethylethylenediamine (TEMED), all stored at 4 °C.

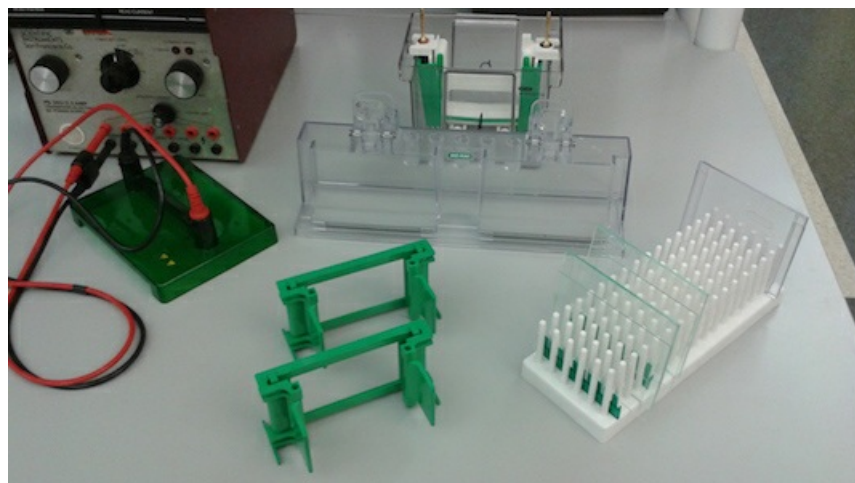

- 2 To assemble the gel casting apparatus, place a short plate on top of a spacer plate. Note that there is an internal space between the two and the writing on the spacer plate is exposed at the top of the sandwich.

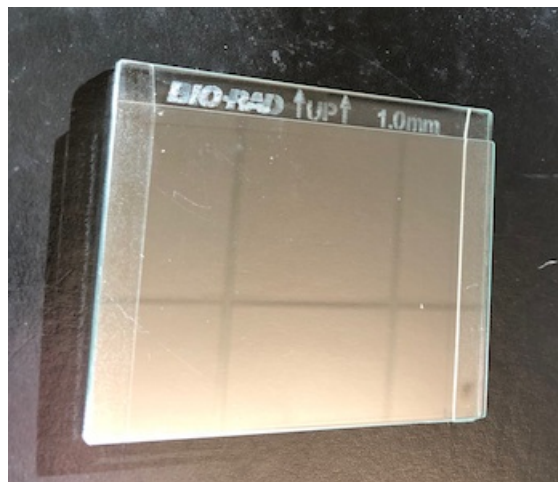

Slide the two plates into an unlocked casting frame (clamps forward) from the top while on a level surface, making sure the short plate is oriented to the front of the frame. With the plates and frame flush with the level surface, lock the assembly by rotating the clamps 90° from their forward to their outward position. Check by hand to ensure both plates and frame are flush at the bottom. Repeat procedure if not.




Place the entire assembly onto a Bio-Rad casting frame with the short plate forward and the casting frame seated with its base back touching the casting stand. Lock the plate/casting stand in place by engaging the spring-loaded lever on the top of the casting stand and applying pressure to the top of the spacer plate with the short plate oriented forward. Now it is ready for gel casting.



- 3 To cast two minigels, add in order: 7.8 mL water, 1.2 mL 5X TBE, 3.0 mL 38:2 acrylamide:bis, and 45 μ L 10% APS to a 15 mL screw cap centrifuge tube. Mix by vortexing briefly. Once everything is ready, add 15 μ L TEMED and vortex briefly to mix before loading gel solution into each of the two casting assemblies. A 6" Pasteur pipette works well for this purpose. Stop before filling (2 mm from top). Carefully insert 10-well comb, using a slight angle to avoid trapping any air, and seat completely. Free-radical polymerization of these gels should be complete in 15 min at room temperature, and these minigels should be ready for use immediately after that. Note: aforementioned 38:2 acrylamide:bis volume is for casting two 10% polyacrylamide gels. Should a different final acrylamide concentration be desired, adjust it and water volumes accordingly.




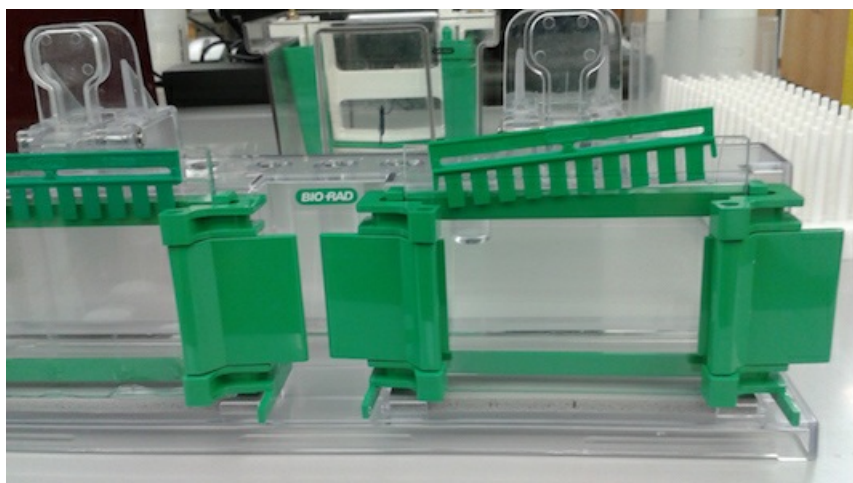


- 4 Note: polyacrylamide minigels in their casting stands can be kept at room temperature/humidity for several hours without any problem. Long-term storage is best when hydrated with a small quantity of water (200 μ L on comb/small plate junction), wrapped in plastic and/or sealed in an airtight container, and stored at 4 °C. Such minigels have effective lifetimes of at least 1 month if stored properly.


Polyacrylamide gel electrophoresis

- 5 With care, secure one polyacrylamide minigel and one buffer dam to each side of the Mini-PROTEAN® electrode assembly. Ensure the green gasket "ears" fully seal with the small plate/spacer junction of the minigel and corresponding plastic lip of the buffer dam before closing side clamps. Place assembly into the tank. Note: we routinely use smaller Mini-PROTEAN® II tanks for buffer economy; adjust subsequent volumes accordingly. Prepare 500 mL $\frac{1}{2}$ X TBE buffer. Fill the central chamber with $\frac{1}{2}$ X TBE buffer and observe if any leakage occurs. If so, drain and repeat seating minigel and buffer dam. Note: if two gels are to be electrophoresed simultaneously, substitute buffer dam for a second minigel.

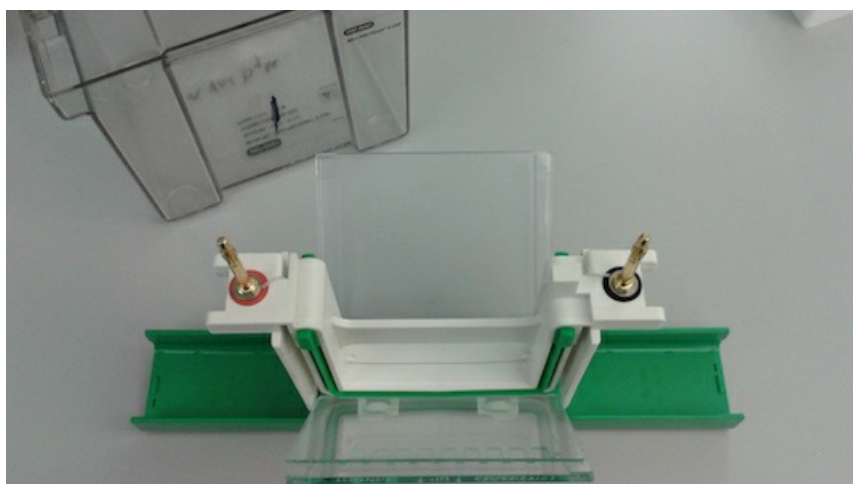


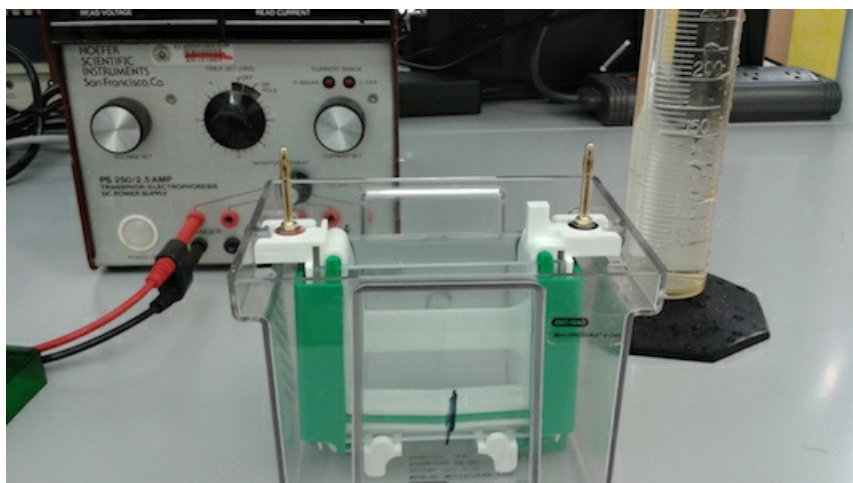

- 6 Prepare DNA samples for electrophoresis. Total volume should not exceed 12 μL and contain at least 3.3% glucose and 0.025% Orange G dye final. **Note:** do not use blue or cyan loading dyes, as many fluoresce in the infrared. Thus, for a 5 μL binding reaction, we add 1 μL 6X Orange Loading Dye (20% glucose, 0.15% Orange G, stored -20°C), and mix thoroughly by repeated pipetting.
- 7 Samples are loaded in order into their respective wells using a P10 Pipetman and 0.1 - 10 μL standard tips. Once all samples are loaded, $\frac{1}{2}\text{X}$ TBE buffer is added to the bottom chamber, the tank lid seated, and leads attached to a power supply. **Note:** be aware of lid-to-electrode assembly and lead-to-power supply socket orientations to maintain proper electrode polarity.

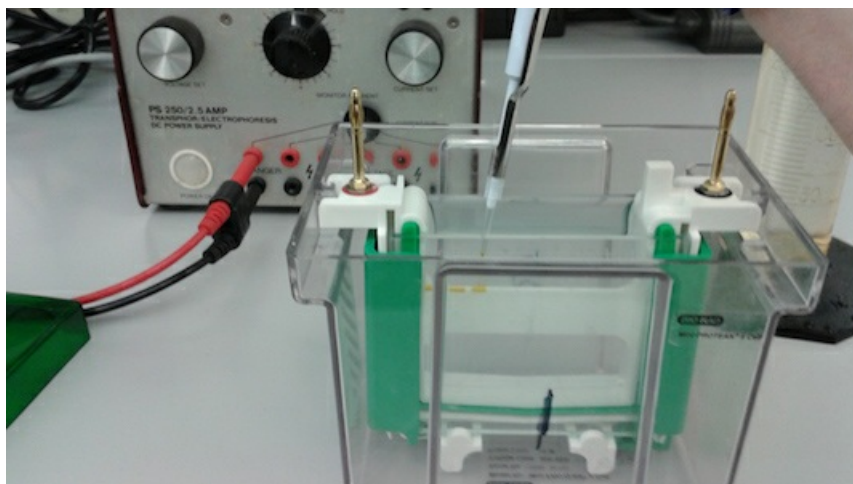


- 8 Electrophorese samples at constant voltage 5 V/cm (50 V) for 5 min to allow the DNA to enter gel, then increase the voltage to 10 V/cm (100 V) for the duration of the run. One hour is sufficient to deliver the Orange G dye to near the bottom of the minigel, which corresponds to approximately 10 bp dsDNA. Once the run is completed, drain tank and electrode assembly, wash with water and disassemble. **Do not take apart gel**, as they can be directly imaged while between glass plates. Note: use of a power supply with integrated timer is a helpful convenience, preventing sample overruns and facilitating reproducibility between runs. However, do not let gels remain at room temperature for extended periods of time, as diffusion of shorter DNAs can occur.

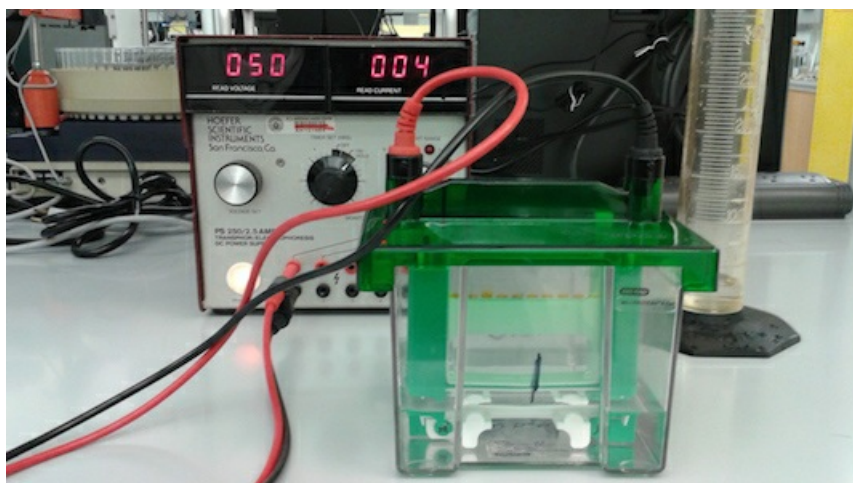


Infrared fluorescence imaging

- 9 To perform infrared fluorescence (IRF) imaging of PAGE minigels, we use a LI-COR Odyssey imager and LI-COR Image Studio™ 5.2 software. In Image Studio™, the DNA preset is used with Scan Controls settings: Scan Resolution (μ): 84 μ m, Scan Quality (Q): Medium, Focus Offset (\oplus): 1.5 mm, Flip Image. A 7 \times 9 cm grid is selected corresponding to the minigel, typically the leftmost, bottom corner of the available field. Check the appropriate channels for the IRDye®-labeled DNAs present. Initial Intensity settings will depend on the quantity of labeled DNA present in each reaction. We use the general guideline: 10 fmoles (Intensity 9), 100 fmoles (Intensity 6), 1000 fmoles (Intensity 3).



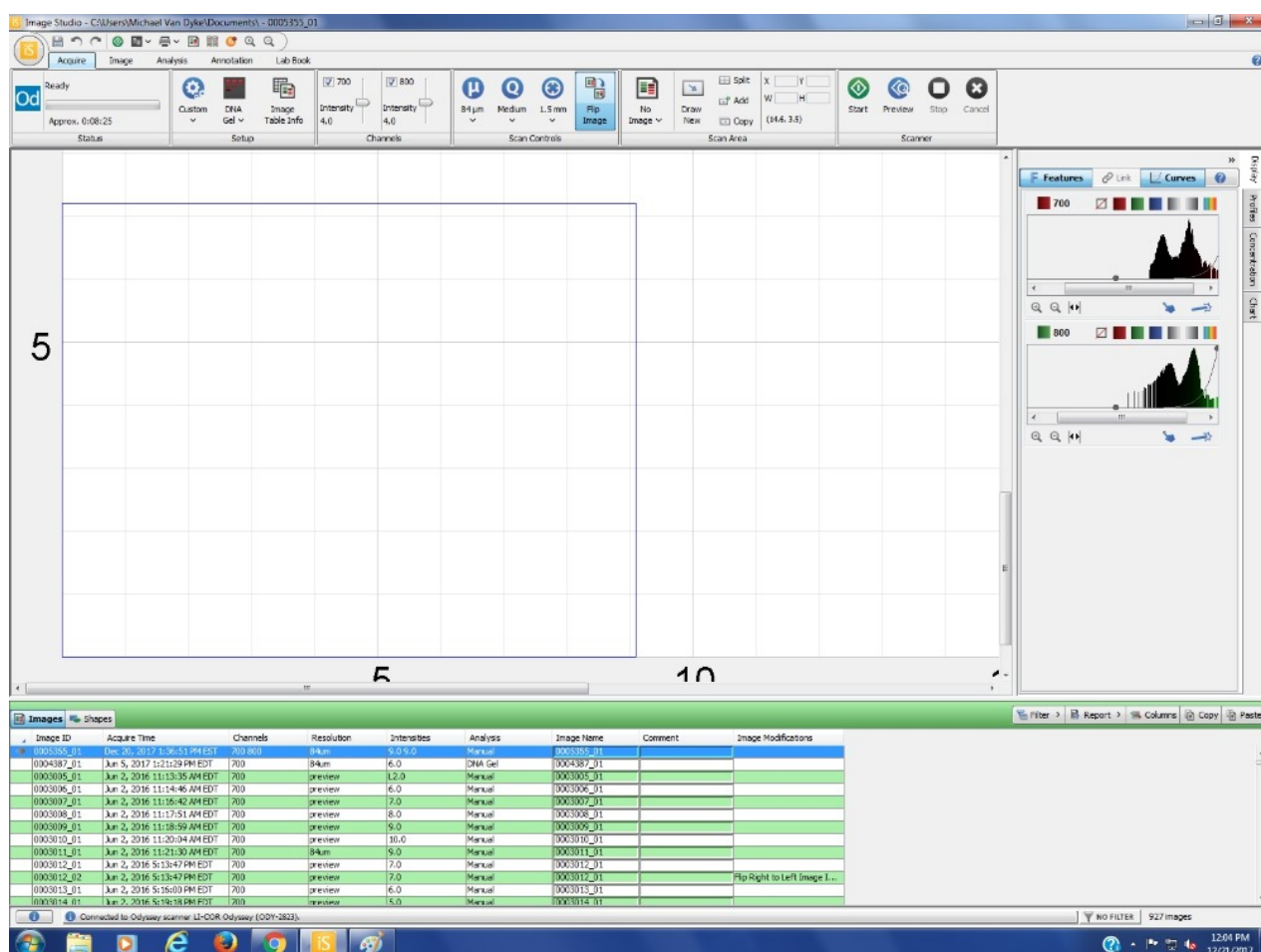


- 10 Note: before imaging, make sure the glass surface of the imager is completely clean. Use DI water and Kimwipes™ to ensure minimal background spotting when should the gel be imaged at high intensity or subjected to high gain enhancement. Place the clean gel sandwich onto the glass imager surface with the short plate face down. Align with both axis and ensure the gel location corresponds to that indicated in Image Studio™. Close imager lid.

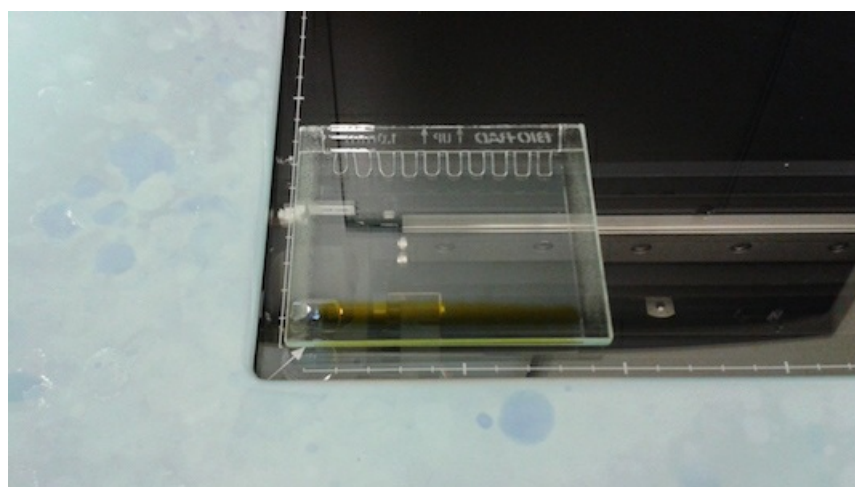



- 11 Initially, a Preview scan at Intensity 4 for both 700 and 800 channels (as needed) is performed, which takes approximately 30 s to run. Additional Preview runs at increasing or decreasing intensity are performed, as needed, until no saturated pixels (white) are observed. Once

maximal Intensity values have been determined, a full run may be initiated (Start). Such takes approximately 8.5 min to run.

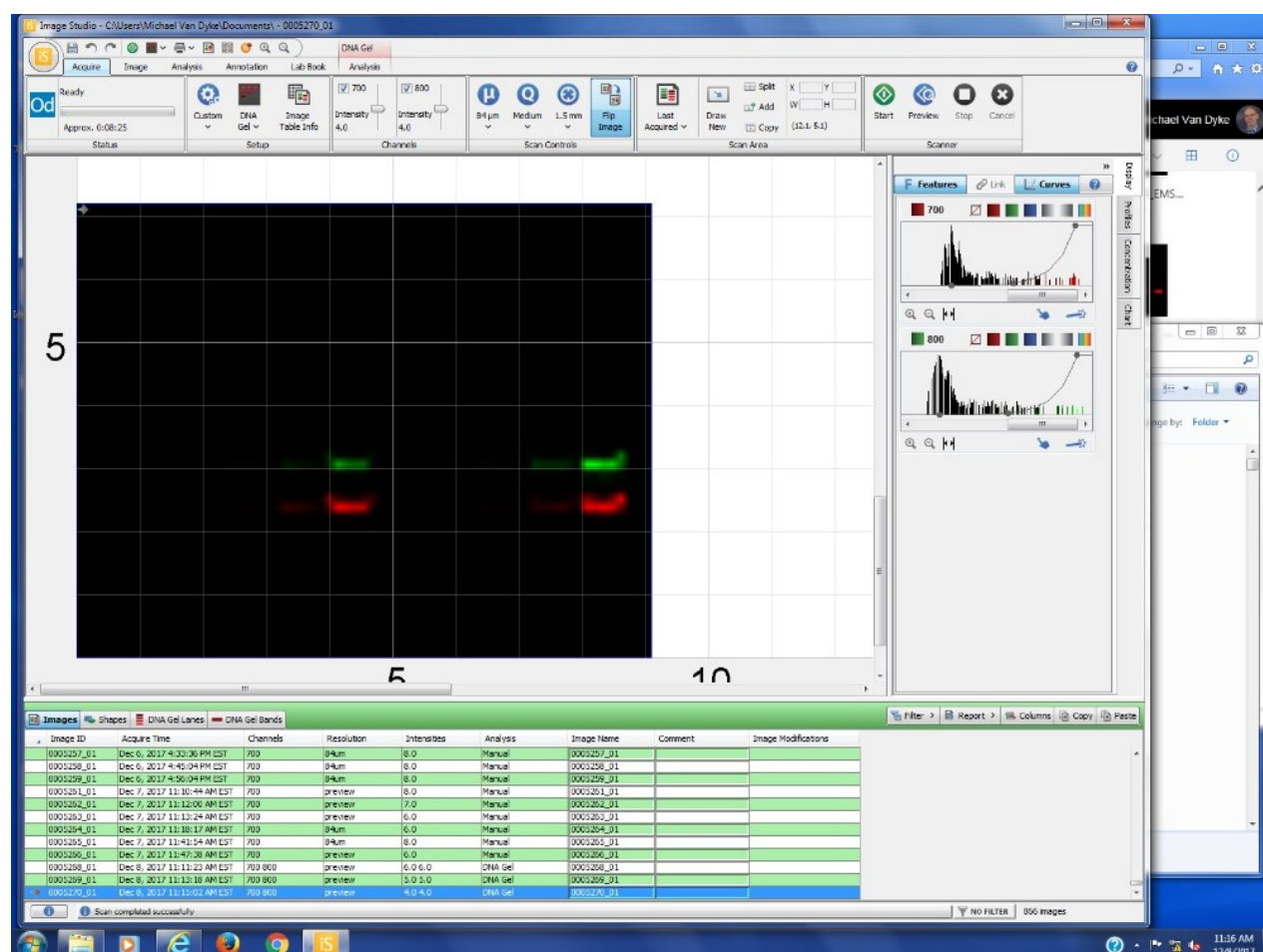


- 12 Images can be saved as tif, png, jpg files and at 150, 300, 600 dpi resolution. Our standard is a 600 dpi jpg file. Gain can be adjusted for each channel using the Curves box, at right. While this allows observation of weaker bands, it is an image manipulation that may not be suitable for publication.

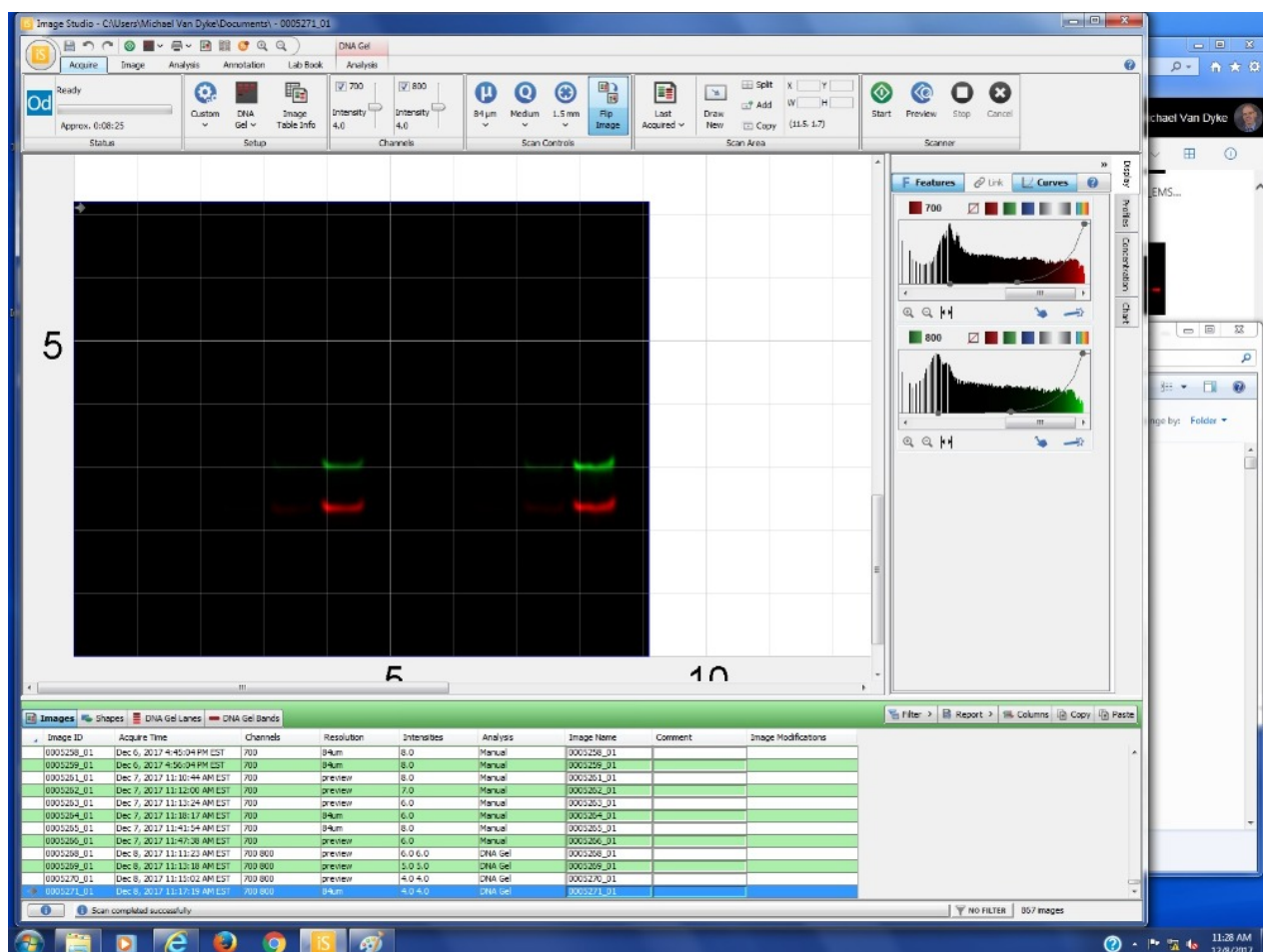


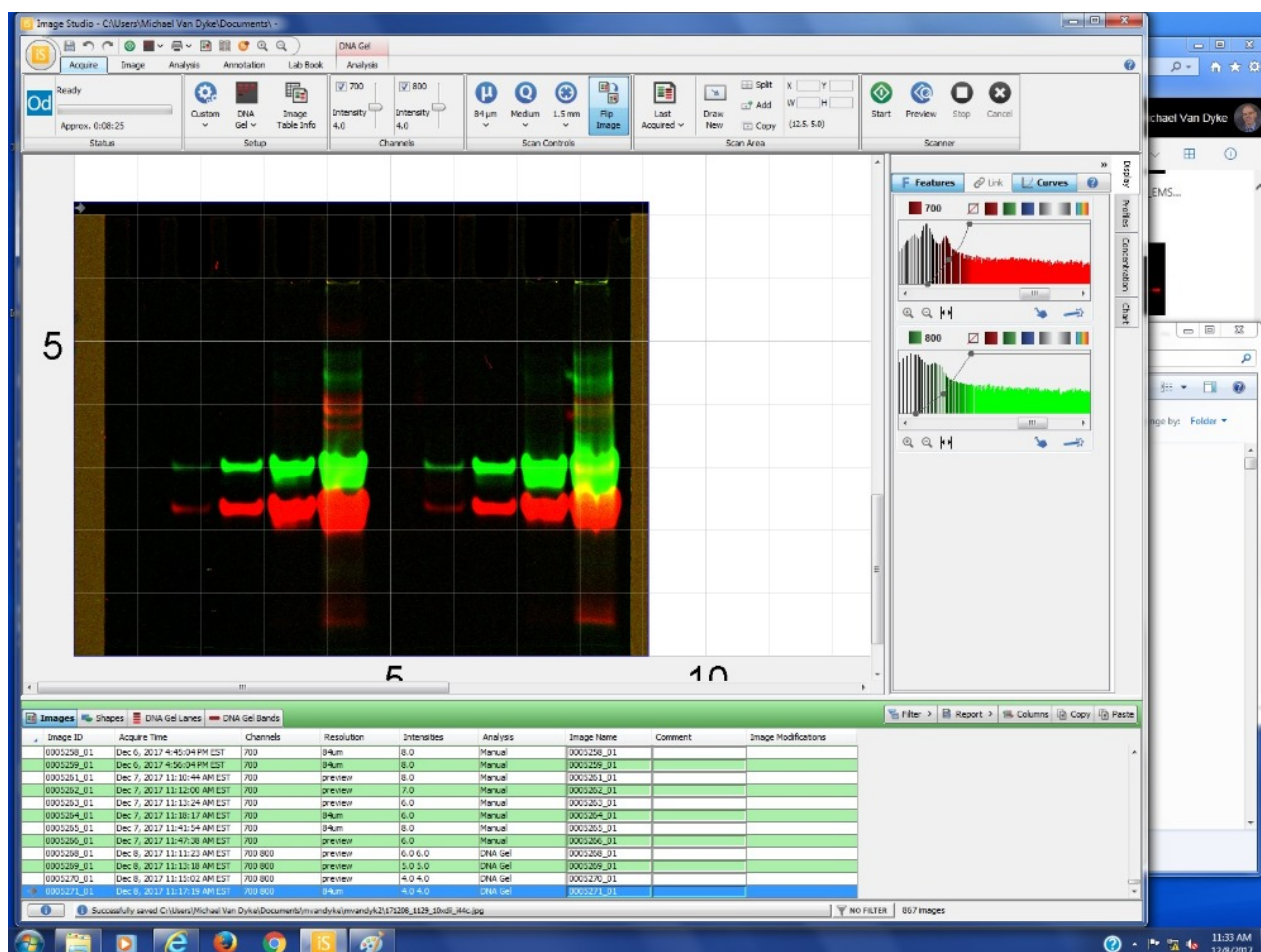


Infrared fluorescence quantitation

- 13 Image Studio™ software provides the capability to quantitate image bands. Note that gain adjustments do not affect these values, so long as no saturated (white) pixels are present. Bands may be automatically identified using the DNA Analysis tab or manually, using the Analysis tab. To use the DNA Gel:Analysis function: (1) Click on the appropriate tab. (2) Click on Lanes:Redraw Boundary to draw a rectangular boundary around the bands of interest. (3) Indicate the number of lanes present. Our default is 10. (4) Click on Bands:Find to find those bands you wish to quantitate. A gray, equivalent sized rectangle will be drawn around each chosen band. Click on Bands:More to identify additional bands for quantitation, as needed.

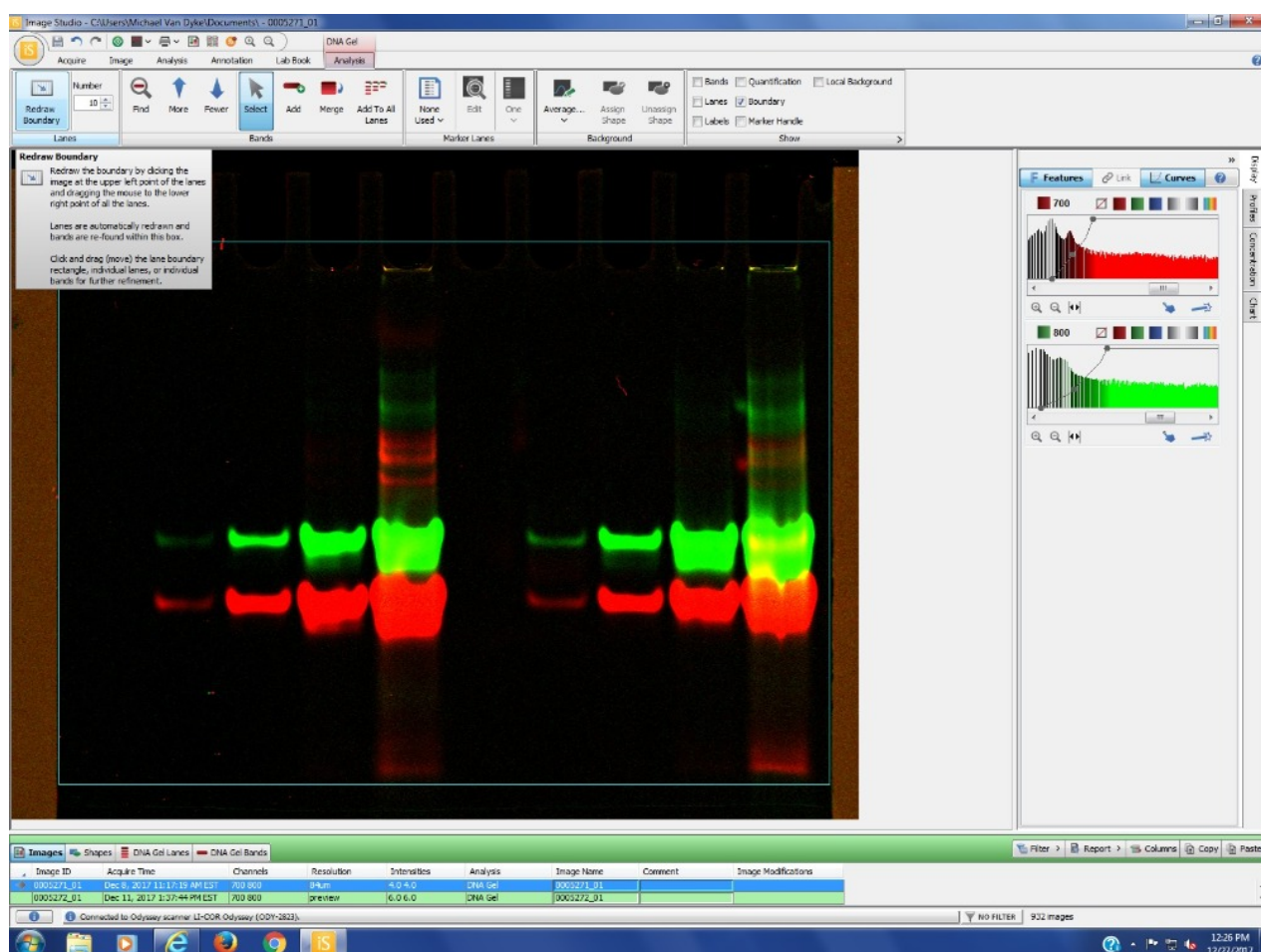


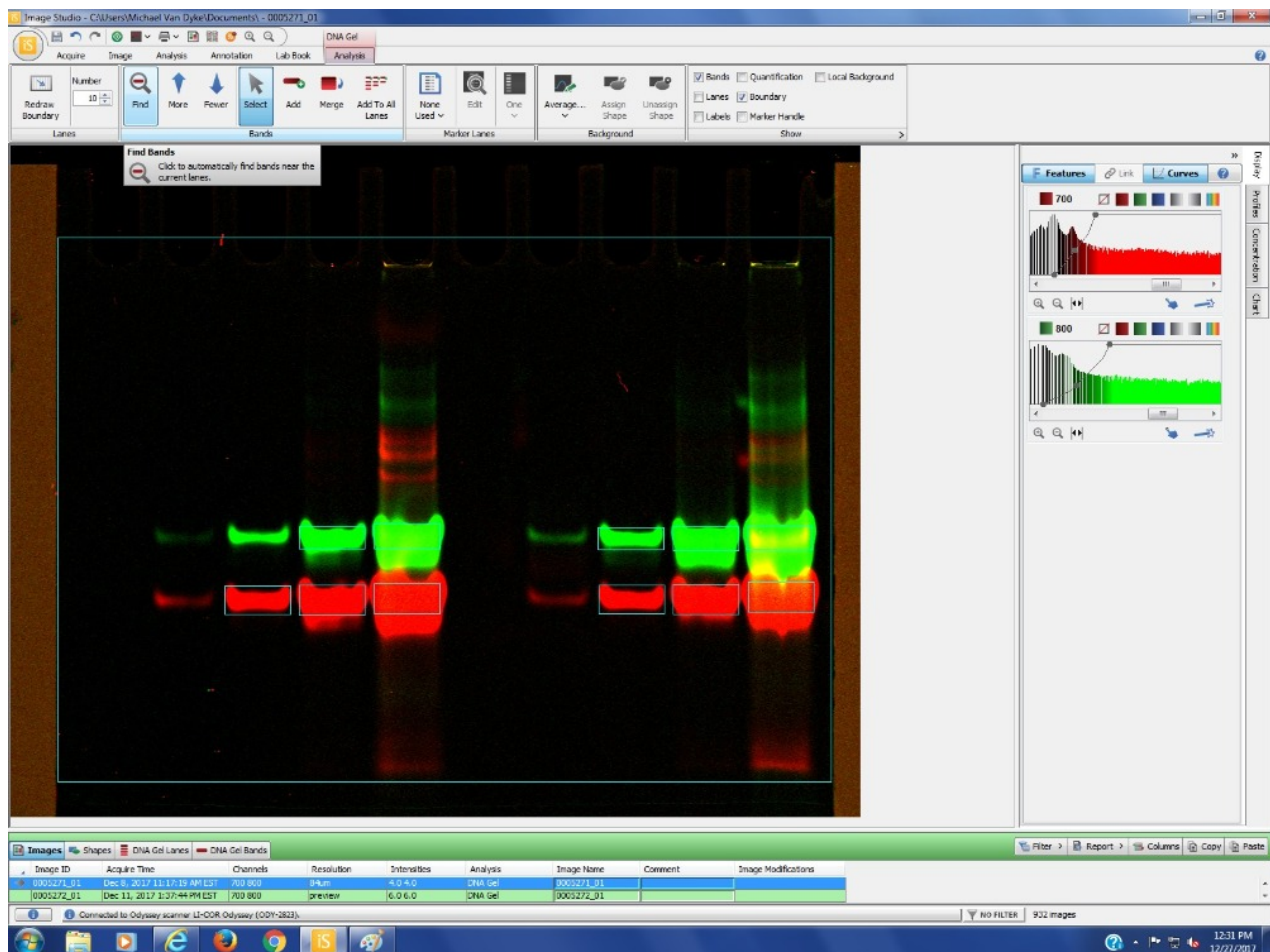

- 14 (5) Once the full complement of bands has been identified, click on Show:Quantification to obtain the pixel numbers for each quantitated band. Note the color bar under each value as these indicate the channel (700 or 800) being quantitated. (6) Remember to save a final image including quantitation values for your records. Occasionally, channel gain (Curves) may need to be reduced to obtain clear images of the pixel numbers. Note that pixel numbers do not change with changes in channel gain.

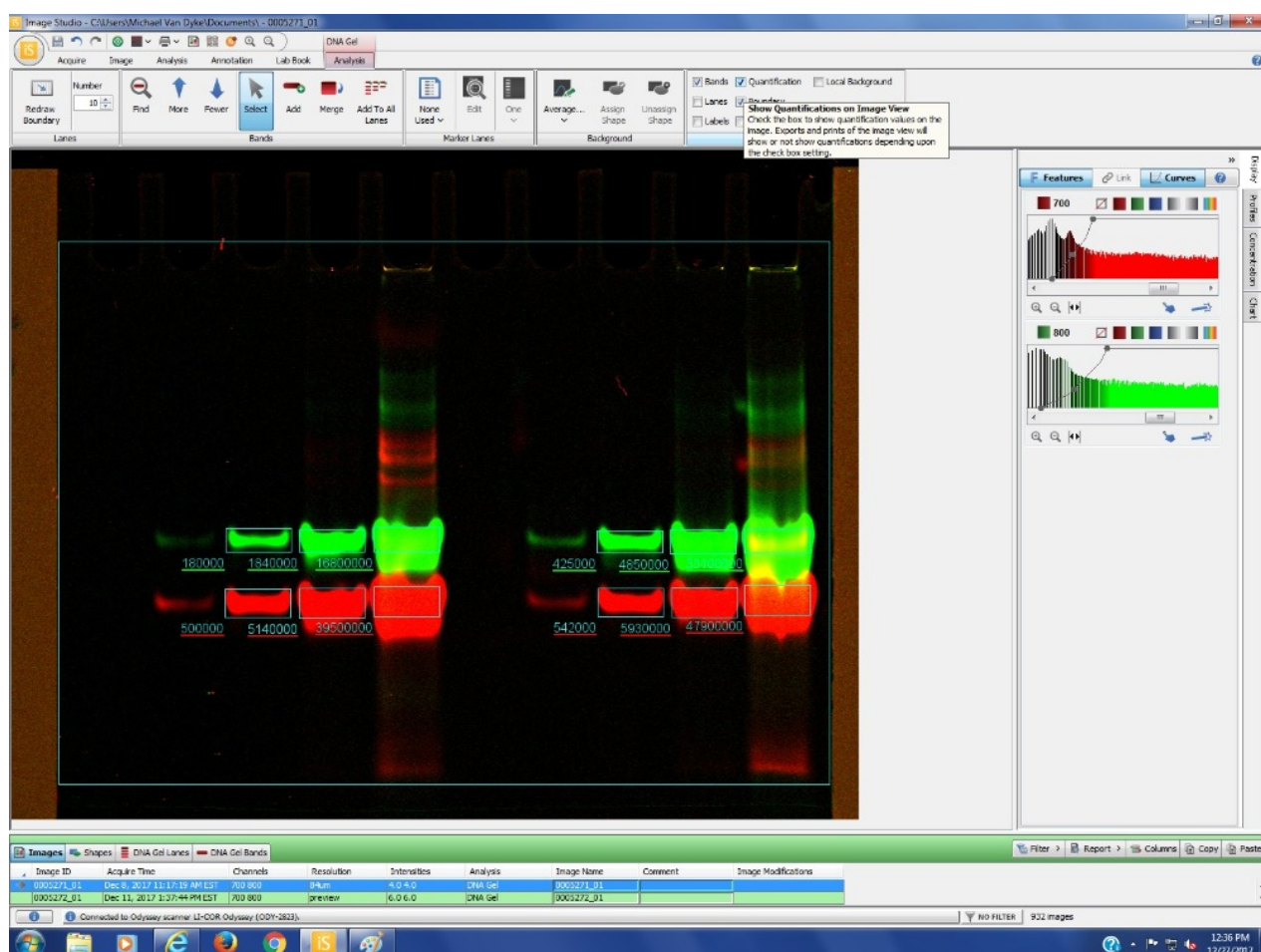

- 15 Under some circumstances (e.g., quantitation of faint bands or background), it is necessary to perform band quantitation manually. Such may be accomplished clicking on the Analysis tab. Use the Shape:Draw Rectangle function to enclose additional bands of interest or background. Take care to ensure all rectangles are of equal dimensions. Note: Analysis requires identifying bands one channel at a time, although both channel analyses may be shown simultaneously once completed. Remember after quantitation to save final images, thereby providing a record of pixel numbers for each band.

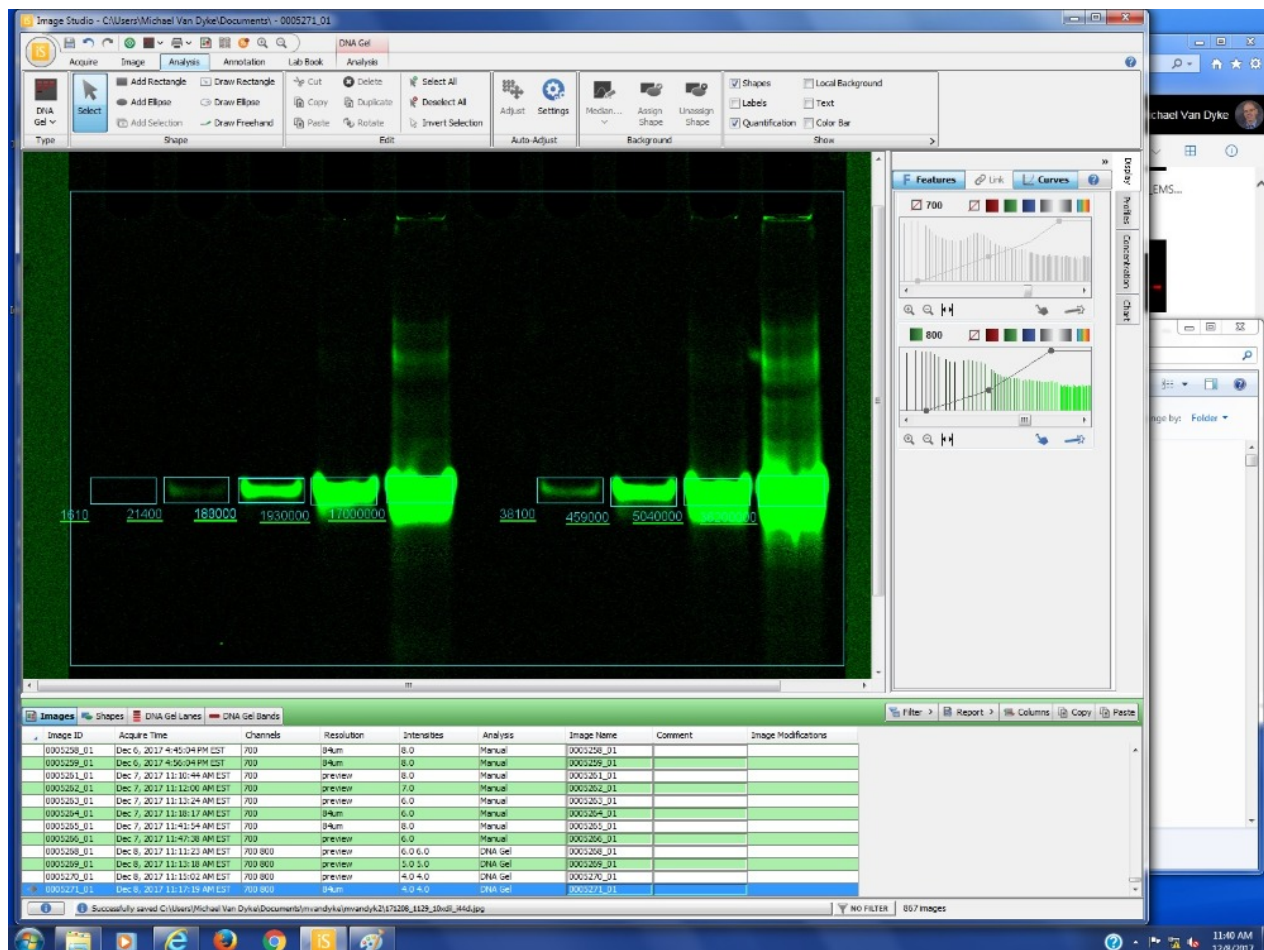

This is an open access protocol distributed under the terms of the [Creative Commons Attribution License](https://creativecommons.org/licenses/by/4.0/), which permits unrestricted use, distribution, and reproduction in any medium, provided the original author and source are credited